

# SPECTROPHOTOMETER CM-3700A

# THE NEW LEADER

Konica Minolta's flagship model: High accuracy and repeatability even at low reflectance/transmittance





# A high-accuracy, top-class model filled with Konica Minolta's advanced optical technology. Ideal for demanding customers with strict requirements for R&D, QC, and CCM applications.

Full data compatibility with CM-3700d

An advanced spectrophotometer for reflectance and transmittance measurements of a broad range of subjects including pigments, dyes, plastic, textiles, paints, ceramics, etc.

# **Top-of-the-line model providing** high accuracy and high reliability

**Konica Minolta's** advanced optical, sensing, and signalprocessing technology provide excellent repeatability.


Strictly selected highquality parts ensure long-term stability and reliability.


**Strict accuracy control** traceable to national standards ensures high quality with high interinstrument agreement.

- 1 Illumination/viewing geometry conforms to CIE, 5 6 Polychromatic unit ISO, ASTM, DIN, and JIS requirements for di:8°, de:8° (diffuse illumination/8° viewing) geometry for reflectance and to CIE, ASTM, DIN, and JIS requirements for transmittance.
- 2 Pulsed xenon arc lamp light source provides high stability, long life, and excellent repeatability even on dark and high-chromaticity colors.
- 3 6-inch integrating sphere has a powdered barium sulfate (BaSO<sub>4</sub>) coating with superior optical characteristics
- 4 Double-beam feedback system directly monitors the light emitted by the xenon lamp at the time of measurement and automatically compensates for changes in brightness or spectral characteristics to ensure high-accuracy measurements.

A diffraction grating S provides high-efficiency separation of light by wavelength, ensuring high repeatability even when measuring dark colors, and the silicon photodiode array sensor 6 quickly converts the separated light into electrical currents. These elements are mounted on stainless steel with a low coefficient of thermal expansion to ensure long-term stability.

- 7 The optical systems for sample measurement and light-source monitoring are mounted on an aluminum alloy block for long-term stability.
  - \* CM-3700A is computer-controlled. Software such as optional SpectraMagic NX2

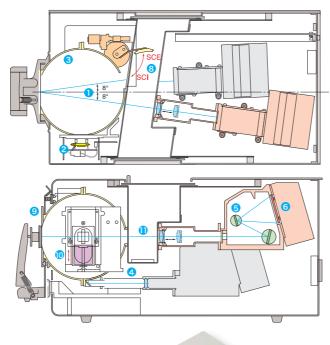

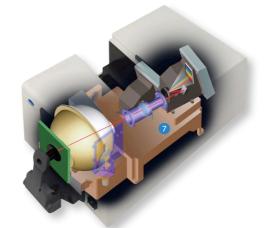

# Top standard instrument for color control systems

With its high accuracy, high repeatability, and high reliability, the CM-3700A can be used as the top standard instrument to which other instruments are referenced in color quality control systems including multiple instruments, both within a company or between companies.



## **Strict quality assurance system** ensures reliability and peace of mind

Konica Minolta's color-measuring instruments are traceable to national standards for wavelength accuracy and white calibration for strict accuracy control. Konica Minolta has also received ISO 9001 and ISO 14001 certification for its integrated quality control system from product development through manufacturing and after-sales service.

### **Comprehensive** support system

Konica Minolta has service facilities worldwide to perform quick inspection, calibration, and repair and ensure that your instruments always provide their best performance.

# **Comprehensive measurement functions**

#### **Reflectance measurements**

8 Switchable between SCI and SCE measurements

SCI (specular component included) measurements minimize the influence of surface conditions on measured values, making it suitable for CCM applications. SCE (specular component excluded) measurements correspond closely to professional visual evaluation

9 Selectable measurement areas

Measurement areas of Ø25.4mm, Ø8mm, and 3×5mm (rectangular) can be selected according to the application.

### 10 Variable UV for measuring fluorescent samples

The UV cut filter can be adjusted in 1000 steps for measurements of samples containing optical whitening agents such as paper, pulp, etc.



#### **Transmittance measurements**

Sideless transmittance chamber for unlimited sample length (Maximum thickness: Approx 50 mm)


di:0°, de:0° (diffuse illumination/0° viewing angle) geometry for spectral transmittance measurements of sheet samples such as glass, filters, etc. as well as cells containing liquid samples such as





#### **Additional Features of CM-3700A**

Long sample holder arm enables measurement at center of A4-size sample.

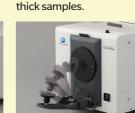


Sample holder arm stays Frame around power open when opened fully switch prevents switching for easy positioning of power off by mistake.



**USB** communication interface for easy connection and highspeed communication



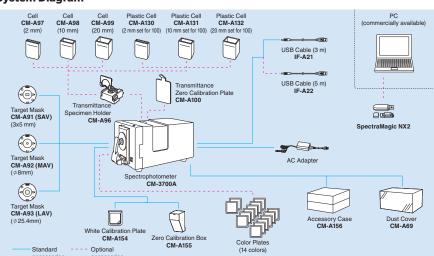

#### **Optional Accessories**

Transmittance Specimen Holder CM-A96



Glass Cell

Transmittance Zero Calibration Plate CM-A100

Color Plates (14 colors)

CM-A97/CM-A98/CM-A99






#### Specifications

Illumination/viewing system	Reflectance	di:8°, de:8° (diffuse illumination/8° viewing angle)				
		SCI (specular component included)/SCE (specular component excluded) switchable				
		Conforms to CIE No. 15 (2004), ISO 7724/1, ASTM E 1164, DIN 5033 Teil 7 and IIS Z 8722 condition c standard.				
	Transmittance					
	- I anomicane	Conforms to CIE No. 15 (2004), ASTM E 1164, DIN 5033 Teil 7 and				
		JIS Z 8722 condition g standard.				
Detector	Silicon photodiode array with flat holographic grating					
Wavelength range	360 to 740 nm					
Wavelength pitch	10 nm					
Half bandwidth	Approx. 14 nm average					
Measuring range	0 to 200%; Resolution: 0.001%					
Light source	Pulsed xenon arc lamp					
Integrating sphere size	ø152 mm					
Minimum measurement interval	3 seconds					
Measurement/illumination area	Reflectance:Changeable between SAV, MAV, and LAV SAV : 3x5 mm measurement / ø5x7 mm illumination MAV : ø8 mm measurement / ø11 mm illumination LAV : ø25.4 mm measurement / ø28 mm illumination Transmittance: Approx. ø20 mm / ø25 mm					
Repeatability	When white calibration plate is measured 30 times at 10-sec. intervals after white calibration has been performed: Spectral reflectance: Standard deviation within 0.05% Chromaticity: Standard deviation within $\Delta E^*$ <sub>ab</sub> 0.005					
	When black tile (BCRA Series II; reflectance: 1%) is measured 30 times at 10-second intervals after white calibration has been performed:  Spectral reflectance: 380 to 740 nm: Standard deviation within 0.02%  360 and 370 nm: Standard deviation within 0.04%  Chromaticity: Standard deviation within △E* <sub>ab</sub> 0.05					
Inter-instrument agreement	Mean $\Delta E^*_{ab}$ 0.08 (typical) Average for 12 BCRA Series II color tiles. Max $\Delta E^*_{ab}$ 0.3 (corresponds to approx. $\Delta E^*_{CMC}$ 0.2) for any of 12 BCRA Series II color tiles compared to values measured with Konica Minolta master body.					
UV adjustment	Computer controlled: continuously variable					
Transmittance chamber	Maximum sample thickness: Approx. 50 mm  Maximum sample length: Unlimited (no sides when transmittance chamber cover is open)  Sample holder (optional) for holding sheet samples or containers of liquid samples can be installed / removed					
	USB 1.1					
Interface						
Interface		/ 50/60 Hz 25 VA (using included AC Adapter)				
Power Operation temperature/	AC 100 to 240 V	<sup>7</sup> 50/60 Hz 25 VA (using included AC Adapter) tive humidity 80% or less (at 33°C) with no condensation				
Power	AC 100 to 240 V 13 to 33°C, rela					
Power Operation temperature/ humidity range (*1) Storage temperature/	AC 100 to 240 V 13 to 33°C, rela 0 to 40°C, relati	tive humidity 80% or less (at 33°C) with no condensation				

#### SpectraMagic NX2 (Optional)

SpectraMagic NX2 is color management software that gives users a customizable screen display and a wide range of functions for operating and configuring their spectrophotometer or Chroma Meter from a computer. Users can display data lists and create color difference graphs and spectral graphs to assist in color management that requires judgment based on numerous values and indicators.

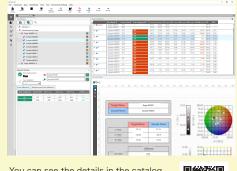

You can see the details in the catalog from the following 2D code. -

SpectraMagic NX2 web Site





- $^{\star}1 \quad \text{Operation temperature/humidity range of products for North America}: 13 \text{ to } 33^{\circ}\text{C}, \text{ relative humidity } 80\% \text{ or less (at } 31^{\circ}\text{C})$ with no condensation
- Displays shown are for illustration purposes only.
- The specifications and appearance shown herein are subject to change without notice.
- without notice. Company names and product names used herein are trademarks or registered trademarks of their respective companies. KONICA MINOLTA, the Konica Minolta logo and symbol mark, "Giving Shape to Ideas", and Spectra Magic are registered trademarks or trademarks of KONICA MINOLTA, INC.

#### SAFETY PRECAUTIONS

For correct use and for your safety, be sure to read the instruction manual before using the instrument.

 Always connect the instrument to the specified power supply voltage. Improper connection may cause a fire or electric shock.



KONICA MINOLTA, INC.	Osaka, Japan					
Konica Minolta Sensing Americas, Inc.	New Jersey, U.S.A.	PHONE: (888)473-2656 (in USA),	+1(201)236-43	00 (outside USA) FAX	: +1(201)785-2	2480 E-Mail: service.sus@konicaminolta.com
Konica Minolta Sensing Europe B.V.	European HQ/ BENELUX German Office French Office UK Office Italian Office Swiss Office Nordic Office Polish Office	Nieuwegein, Netherlands München, Germany Roissy CDG Cedex, France Warrington, United Kingdom Cinisello Balsamo, Italy Dietikon, Switzerland VÄSTRA FRÖLUNDA, Sweden Wrocław, Poland	PHONE: PHONE: PHONE: PHONE: PHONE: PHONE: PHONE: PHONE:	+31(0)30 248-1193 +49(0)89 4357 1560 +33(0)1 80 11 10 70 +44(0)1925 467300 +39 02849488.00 +41(0)43 322-9800 +46(0)31 7099464 +48(0)71 73452-11	E-Mail: E-Mail: E-Mail: E-Mail: E-Mail: E-Mail: E-Mail: E-Mail:	info.benelux@seu.konicaminolta.eu info.germany@seu.konicaminolta.eu info.france@seu.konicaminolta.eu info.tuk@seu.konicaminolta.eu info.uk@seu.konicaminolta.eu info.sitaly@seu.konicaminolta.eu info.sitaly@seu.konicaminolta.eu info.mordic@seu.konicaminolta.eu info.poland@seu.konicaminolta.eu info.poland@seu.konicaminolta.eu
Konica Minolta (CHINA) Investment Ltd.	SE Sales Division Beijing Office Guangzhou Office Chongqing Office Qingdao Office Wuhan Office Shenzhen Office Xi'an Office	Shanghai, China Beijing, China Guangzhou, China Chongqing, China Shandong, China Hubei, China Shenzhen, China Xi'an, China Xiamen, China	PHONE: PHONE: PHONE: PHONE: PHONE: PHONE: PHONE:	+86-(0)21-6057-1089 +86-(0)10-8522 1551 +86-(0)20-3826 4220 +86-(0)23-6773 4988 +86-(0)532-8079 187 +86-(0)27-6885 0586 +86-(0)755-2868 753 +86-(0)592-7107 399	E-Mail: E-Mail: E-Mail: 1 E-Mail: E-Mail: 5 E-Mail: E-Mail:	hcn_sensing@gcp.konicaminolta.com hcn_sensing@gcp.konicaminolta.com hcn_sensing@gcp.konicaminolta.com hcn_sensing@gcp.konicaminolta.com hcn_sensing@gcp.konicaminolta.com hcn_sensing@gcp.konicaminolta.com hcn_sensing@gcp.konicaminolta.com hcn_sensing@gcp.konicaminolta.com
Konica Minolta Sensing Singapore Pte. Ltd.	Singapore		PHONE:	+65 6563-5533	E-Mail:	se-service.sg@konicaminolta.com
Konica Minolta Sensing Korea Co., Ltd.	Korean HQ Cheonan Office	Goyang-si, Korea Cheonan-si, Korea	PHONE: PHONE:	+82(0)2-523-9726 +82(0)41-556-9726	E-Mail: E-Mail:	se.korea@konicaminolta.com se.korea@konicaminolta.com