

# **Spectrophotometer**

## **CM-26dG**

## **CM-26d**

## **CM-25d**

### **En Basic Operating Instructions**



Please read before using the instrument.

This manual explains the basic procedures for measuring color and/or gloss with the CM-26dG / 26d / 25d spectrophotometers. Use this 2D code to download this manual in pdf form.

Note)

For setting and operation details, see the instruction manual of the CM-26dG / 26d / 25d.



KONICA MINOLTA


# Table of Contents

Names of Parts .....	1
Measurement .....	5
When color difference is measured .....	9
Setting measurement conditions .....	16
How to set the measuring instrument .....	17

# Names of Parts

## ■ Instrument

**Viewfinder  
(Specimen  
confirmation  
window)**

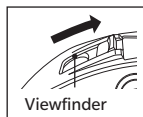

**Viewfinder lever**

**LCD screen**

**Control panel**

**Charging lamp**

**Power switch**


**USB connection  
terminal  
(Mini-B type)**

**Measurement button  
(It is on the left and right)**

**Wrist strap  
attachment hole**

**Sample  
surface  
aperture**

**Measurement  
area switch**

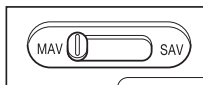


**(CM-26dG/26d only)**

**Battery  
cover**

**Screw holes for securing**


## ■ Display (LCD Screen) P.22, Instruction Manual

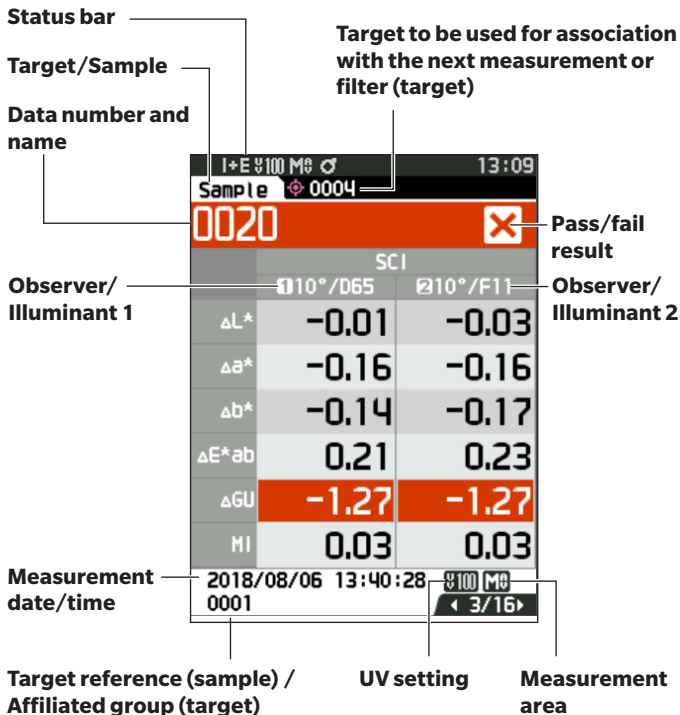

### ■ Operation Keys

**[/↗] (Target/Sample) key**  
Switches between **<Target> screen** and **<Sample> screen**.

**[MENU] key**  
Displays the **<Settings> screen**.

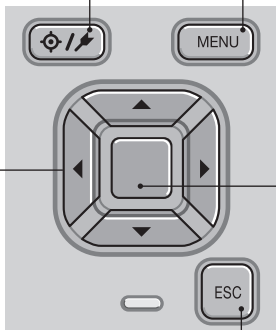


**[Confirmation] key**  
Sets the item or setting indicated by the cursor on the various setting screens. This key also switches to detail screens for data selected in the list displayed on the **<Results Display> screen**.

**[◀, ▶, ▲, ▼] keys**  
Switches screen tabs from the **<Results Display> screen**, moves the cursor on the **<Settings> screen**, or changes the selected value.


**[ESC] key**  
Returns to the previous screen without configuring the settings when pressed on the **<Settings> screen**, and returns to list screen when pressed on the sample details screen.

## ■ Calibration Stage


CM-26dG

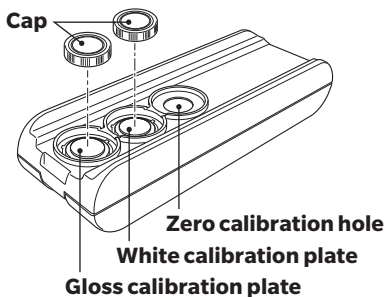

CM-26d

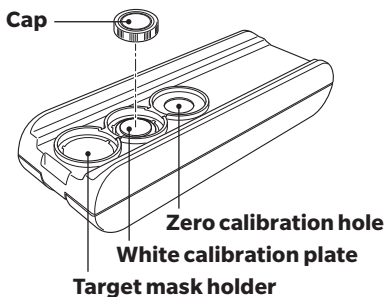

CM-25d

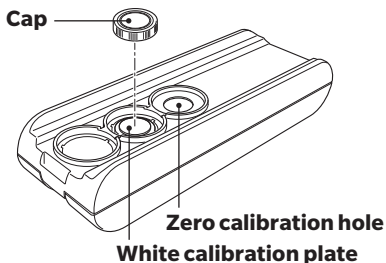

# Measurement

## 1 Turning the Power ON (P.6)



Language selection

**P.110, Instruction Manual**

\* As necessary (after initialization, etc.)



Measurement cond. setting **P.99, Instruction Manual**



Measurement area selection \* When required depending on the measurement conditions

**P.33, Instruction Manual**

**Note: Make sure the measurement area set on the instrument is the same as the area of the target mask to be used.**



Zero calibration

**P.35, Instruction Manual**

\* Only when required (when changing the measurement conditions, etc.)



## 2 White calibration (P.6, P.7)

**White calibration**

**Gloss calibration (\* CM-26dG only)**



## 3 Measurement (P.8)

# 1

## Turning the Power ON

P.32,  
Instruction Manual

- 1 Press and hold the power switch for about 1 second.



- When first turning the instrument ON after purchasing, the language setting screen will be displayed, followed by the date and time setting screen. Configure the settings according to P.103 and P.104.

# 2


## White calibration

P.37,  
Instruction Manual


### To perform white calibration

**Note:** Zero calibration must be performed before white calibration if there have been significant changes in the measurement environment, if the instrument has not been used for a long period of time or if using the MAV target mask (with glass).

- 1 Select "Calibration (Excluding Zero Cal.)" and press the [Confirmation] key.

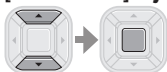



- 2 Place the instrument on the calibration stage.

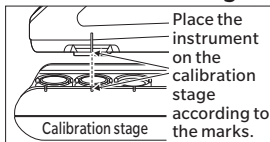

- 3 Press the measurement button.



- White calibration will be performed.
- Do not move the instrument until white calibration is complete.

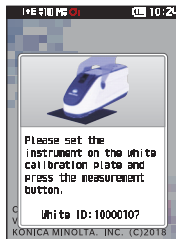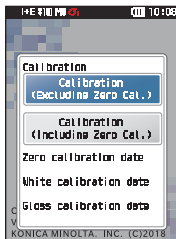


## Gloss calibration only with the 26dG

P.38,  
Instruction Manual

### If using the CM-26dG, perform gloss calibration next.



- ④ **Place the instrument in the gloss calibration position.**

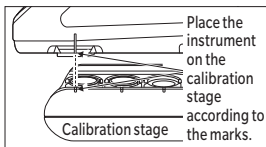

- ⑤ **Press the measurement button.**



- Gloss calibration will be performed.
- Do not move the instrument until gloss calibration is complete.
- When gloss calibration is completed, the screen returns to the <Sample> screen.

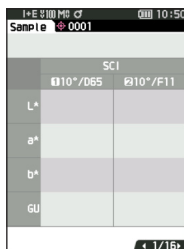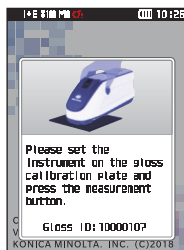


## 3

### Measurement

P.41,  
Instruction Manual

#### ① Check the Sample screen.

- If not displayed, switch to the Sample screen.

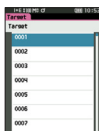

<Reference  
color screen>




<measurement  
screen>

I+E 5100 M0 C 10:54		
Sample 0001		
SCI		
	10°/D65	2°/F11
L*		
a*		
b*		
GU		
1/16		

#### ② Set the specimen measuring port on the sample to measure.

- Take care to prevent floating or tilting from occurring.

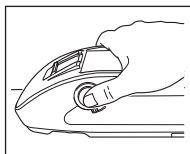

#### ③ Press the measurement button.



- The results are displayed on the screen. The sample was saved as "0001".

I+E 5100 M0 C 10:56		
Sample 0001		
0001		
SCI		
	10°/D65	2°/F11
L*	47.62	52.40
a*	52.42	52.37
b*	27.59	35.57
GU	12.72	12.72
2019/02/20 10:56:19 100 Ms		
0001 1/16		

# When color difference is measured

**<Measuring color difference in the normal mode>**

**1 Color difference target settings (P.10)**



**2 Pass/fail judgment criteria settings (P.11)**



**3 Measurement (P.13)**

**<Measuring color difference in the simple mode>**

**1 Set to "Simple mode". (P.14)**



**2 Measuring a target (P.15)**



**3 Measuring a sample (P.15)**

## <Measuring color difference in the normal mode>

# 1

## Color difference target settings


P.41,  
Instruction Manual

### Setting color difference by measuring a target

#### ① Display the target or target list screen.



- Set the target number with the [▲] and [▼] keys before imparting measurement.

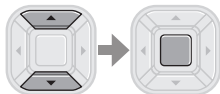


SCI		
	D10°/D65	D10°/F11
L*		
a*		
b*		
GU		

1 / 2

#### ② Set the specimen measuring port on the sample to measure.

- Take care to prevent floating or tilting from occurring.

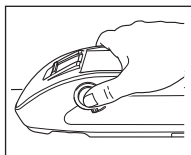

#### ③ Press the measurement button.



- The results are displayed on the screen.
- The selected number is used as the target data number. If data already exists under that number, the instrument asks whether to overwrite that data or not, therefore enter the target by pressing the [Confirmation] key.

SCI		
	D10°/D65	D10°/F11
L*	47.60	52.37
a*	52.52	52.40
b*	26.88	34.79
GU	15.78	15.78

2019/02/20 10:57:39 1 / 2


The target is set.

# 2



## Pass/fail judgment criteria settings

P.60-61, Instruction Manual

- ① Select <Target Menu> - <Pass/Fail> - "Edit tolerance".


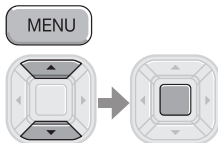


- ② Input settings.

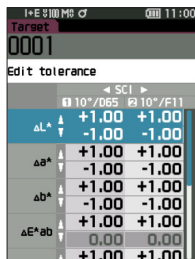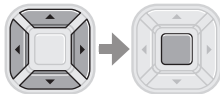


### Settings

- ☐ x / y : -0.2000 to 0.2000
- ☐ Equation / MI : 0.00 to 20.00
- ☐ Other than the above : -20.00 to 20.00

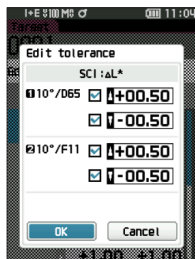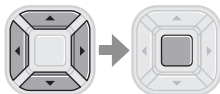

- ③ Once settings have been completed, select "OK" to establish them.

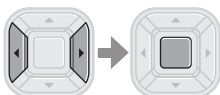

I+E 9100 M0 11:04

Target 0001

Edit tolerance

	SCI	
	10°/D65	10°/F11
$\Delta L^*$	+0.50	+0.50
	-0.50	-0.50
$\Delta a^*$	+1.00	+1.00
	-1.00	-1.00
$\Delta b^*$	+1.00	+1.00
	-1.00	-1.00
$\Delta E^*ab$	+1.00	+1.00
	0.00	0.00
	+1.00	+1.00

- ④ Close the "Edit tolerance" screen.



I+E 9100 M0 11:05

Target

Target menu[0001]

Pass/Fail

Edit tolerance

Tolerance list

Warning level 80%

Parametric coef.

# 3

## Measurement

P.41,  
Instruction Manual

### ① Display the Sample screen.




- Select a target number before imparting measurement.

Sample 0001		11:06
0001		
SCI		
10°/D65		10°/F11
L*	47.62	52.40
a*	52.42	52.37
b*	27.59	35.57
GU	12.72	12.72
2019/02/20 10:56:19		0001

### ② Set the specimen measuring port on the sample to measure.

- Take care to prevent floating or tilting from occurring.

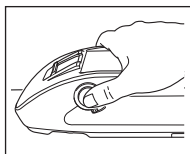

### ③ Press the measurement button.



- The results are displayed on the screen.
- A sample number is automatically assigned in the order of measurement.

Sample 0001		11:08
0002		
SCI		
10°/D65		10°/F11
L*	47.69	52.47
a*	52.44	52.40
b*	27.67	35.62
GU	12.93	12.93
2019/02/20 11:08:17		0001

✓ Mark signifying a "Pass" judgment


# <Measuring color difference in the simple mode>

# 1



## Set to "Simple mode".

P.48,  
Instruction Manual

- ① Press the [ESC] button to skip the calibration prompt displayed immediately after turning the power ON.


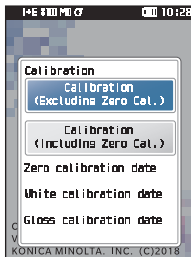


- ② Select "Yes" when "Change to the simple mode?" appears on the display.

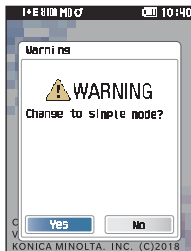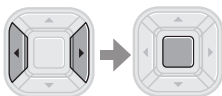

Or,

- ① Select <Settings> - <Instrument Mode> - "Simple".

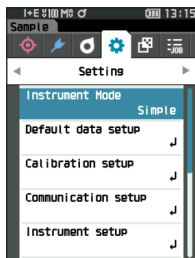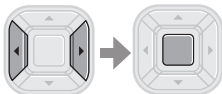

- ② Press the [Target/Sample] button or [ESC] button to switch to the simple measurement screen.



## 2

### Measuring a target:

Absolute value measurement

P.48,  
Instruction Manual

- ① Press the measurement button.



Target		
SCI		
	10°/D65	2°/F11
L*	47.60	52.37
a*	52.52	52.40
b*	26.88	34.79
GU	15.78	15.78
2019/02/20 13:20:39		
1/2		

## 3

### Measuring a sample:

Difference measurement

P.48,  
Instruction Manual

- ① Press the measurement button.



#### Notes

- Measurement data is not saved in the simple mode.
- The opacity mode cannot be selected while the instrument is in the simple mode. Moreover, the simple mode cannot be engaged while the opacity mode is active.

Sample		
SCI		
	10°/D65	2°/F11
ΔL*	0.17	0.16
Δa*	-0.13	-0.05
Δb*	0.43	0.45
ΔE*ab	0.48	0.48
ΔGU	-0.14	-0.14
HI	0.08	0.08
2019/02/20 13:20:59		
3/5		

# Setting measurement conditions

Measurement conditions (Measurement mode, SMT Times, Observer/Illuminant, Display) must be set before imparting measurement.

The menu items that can be set are as follows.

## **P.87 - 107, Instruction Manual**

### Measurement cond.

#### Measurement mode

Color & Gloss / Color only / Gloss only / Opacity

#### Measurement setup

Specular component

UV

Auto average

Manual average

Manual average option (SMC average option)

SMC

SMC Threshold

SMC Times

#### Observer/Illuminant

Observer/Illuminant 1

Observer/Illuminant 2

#### Display cond.

Display type

Color space

Color difference equation

Custom 01 to 14




# How to set the measuring instrument

To set measurement instrument options, select “Instrument setup” from the <Setting> screen.

The menu items that can be set are as follows.

**P.108 - 117, Instruction Manual**

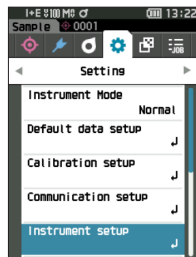

## Instrument setup

User type

Language

Date format

Date & time

Brightness

Direction

Beep

Auto power off

Password setting



KONICA MINOLTA

En 9222-AC5J-43  
CBHAKK