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Installing the software

When installing the software, you must log on the computer as a user with Administrator privileges. If you are using a previous version of Configuration Tool CM-CT1, please uninstall the old Configuration Tool CM-CT1 before installing the new one.

1. Installation procedure

[Windows 10]

1. Double-click the [setup.exe] .
2. Click "Next".
3. Read the software license agreement carefully. Select "I accept the terms in the license agreement" and click "Next".
4. The set-up program asks you to choose the folder in which you want to install the software. To install the software in another location, click the "Change" button, type the path to the location and click "OK".
5. Click "Next" .
6. Click "Install" .

The set-up program will begin copying the files.

7. If dialog for User Account Control appears, click "Yes".
8. This completes the installation process. Click "Complete".

A shortcut is created on the desktop.

[Windows 8.1]

If .NET Frameworks 4.5.2 is installed on your computer, it will be the same as for Windows 10.

1. Double-click the [setup.exe] .
2. As the dialog for starting installation of .NET Frameworks 4.5.2 appears, click "OK" .
3. If dialog for User Account Control appears, click "Yes".
4. After installing .NET Frameworks 4.5.2, the computer will automatically restart.
5. After restarting the computer, installation of this software starts automatically.
6. Click "Next".
7. Read the software license agreement carefully. Select "I accept the terms in the license agreement" and click "Next".
8. The set-up program asks you to choose the folder in which you want to install the software. To install the software in another location, click the "Change" button, type the path to the location and click "OK".
9. Click "Next".
10. Click "Install".

The set-up program will begin copying the files.

11. If dialog for User Account Control appears, click "Yes"
12. This completes the installation process. Click "Complete".

A shortcut is created on the desktop.

2. To uninstall

[Windows 10]

1. Right-click in the lower-left corner of the screen.
2. Click [Application and Features] on the menu that appears.
3. Select [Configuration Tool CM-CT1] , and then click “Uninstall”.
4. A dialog box for confirming that you want to uninstall the application appears. Click “Yes”.
5. A dialog box for confirming that you want to update the software appears. Follow the instructions on the screen to complete the uninstall process.

[Windows 8.1]

1. Right-click in the lower-left corner of the screen.
2. Click [Programs and Features] on the menu that appears.
3. Select [Configuration Tool CM-CT1] , and then click “Uninstall”.
4. A dialog box for confirming that you want to uninstall the application appears. Click “Yes”.
5. A dialog box for confirming that you want to update the software appears. Follow the instructions on the screen to complete the uninstall process.