# Data Management Software

## CS-S20

- E** Installation Guide
- D** Installationsanleitung
- F** Guide d'installation
- I** Guida all'installazione
- Es** Guía de instalación
- CS** 安装指南
- CT** 安装指南
- J** インストールガイド

KONICA MINOLTA

English

Deutsch

Français

Italiano

Español

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## Software License Agreement

The terms of the license agreement of the CS-S20 software are provided in the Software License Agreement dialog box displayed on-screen during the installation process. This software can be installed only if you agree to all the terms of the agreement.

### Safety Precautions

Before you use the CS-S20 software, we recommend that you thoroughly read this manual as well as the instruction manuals of your PC and the instrument.

## Notes on Installation

- The CS-S20 application software is designed to be used with the Windows 7, Windows 8, Windows 8.1 or Windows 10 operating system. Note that neither operating system is included with this software.
- One of these operating systems must be installed on the PC before this software can be installed.
- When inserting the CD-ROM into the CD-ROM drive, note the correct orientation of the disc. Insert it gently.
- Keep the CD-ROM clean and free from scratches. If the recorded surface becomes dirty or the label surface is scratched, a read error may result.
- Avoid exposing the CD-ROM to rapid temperature changes and condensation. Avoid leaving it in locations where it may be exposed to high temperatures from direct sunlight or heaters.
- Do not drop the CD-ROM or subject it to strong impact.
- Keep the CD-ROM away from water, alcohol, paint thinners, and other such substances.
- Remove the CD-ROM from the CD-ROM drive while the computer is turned on.

## Notes on Storage

- After using the CD-ROM, return it to its case and store in a safe place.
- Avoid leaving the CD-ROM in locations where it may be exposed to high temperatures from direct sunlight or heaters.
- The CD-ROM should not be kept in areas of high humidity.
- Every effort has been made to ensure the accuracy of the content of this software. However, should you have any questions or comments, please contact the nearest KONICA MINOLTA authorized service facility.

## Disposal Method

- Make sure that this product (CD-ROM and protection key) is either disposed of or recycled correctly in accordance with local laws and regulations.

## Formal designations of application software used in this manual

(Designation in this manual)	(Formal designation)
Windows, Windows 7	Microsoft® Windows® 7 Professional Operating System
Windows, Windows 8	Microsoft® Windows® 8 Pro Operating System
Windows, Windows 8.1	Microsoft® Windows® 8.1 Pro Operating System
Windows, Windows 10	Microsoft® Windows® 10 Pro Operating System

## Trademarks

- “Microsoft”, “Windows”, “Windows 7”, “Windows 8”, “Windows 8.1”, and “Windows 10” are registered trademarks of Microsoft Corporation in the U.S.A. and other countries.
- Other company names and product names mentioned in this manual are the registered trademarks or trademarks of their respective companies.

## Package Contents

- Installation CD-ROM of CS-S20 (Qty: 1)
- Installation Guide (this manual)
- Authorized Service Facility

The Instruction Manual is also installed in PDF form with a shortcut in the start menu during the software installation.

To read the manual, go to Start Menu → All programs → KONICA MINOLTA → CS-S20 → CS-S20 Manual.

\* For Windows 8 or Windows 8.1, a shortcut to the manual is registered to a different screen. To display this screen, right-click in the Start screen, and then click All apps in the lower right of the screen.

Versions of the Instruction Manual in other languages are also included on the installation CD-ROM.

You will need Adobe Reader<sup>®</sup> from Adobe Corporation. The latest Adobe Reader<sup>®</sup> can be downloaded for free from the Adobe website. Also, it is possible to use the Adobe Reader<sup>®</sup> installer included on the installation CD-ROM.

(Example) When the CD-ROM is in E drive E:\Adobe Reader\EN

# 1. Installing the software

The CS-S20 software is provided on a CD-ROM. A CD-ROM drive is required to install the software. When installing the software, you must log on to the computer as a user with Administrator privileges.

## 1.1 Installation procedure

Install the CS-S20 data management software.

Instrument USB drivers are automatically installed while installing CS-S20.

A dialog box for checking the User Account Control may be displayed multiple times during the process as the software and various drivers are installed. When this dialog box is displayed, click “Continue”.

### [Windows 7]

1. Start Windows on your computer.
2. Insert the CS-S20 CD-ROM into the CD-ROM drive.
3. The install program will start automatically. If it does not start automatically, double-click the [setup.exe] icon. → Proceed to step 4.

### [Windows 8, Windows 8.1 and Windows 10]

1. Start Windows on your computer.
2. Insert the CS-S20 CD-ROM into the CD-ROM drive.
3. On the screen, the message “Tap to choose what happens with this disk.” is displayed. Click this message, and then click [Run setup.exe].
4. The Select Language dialog box opens. Select the appropriate language and click “OK”.  
The following procedure assumes that English is selected.
5. Click “Next”.
6. Read the software license agreement carefully. Select “I accept the terms in the license agreement” and click “Next”.
7. The set-up program asks you to choose the folder in which you want to install the software. To install the software in another location, click the “Change” button, type the path to the location and click “OK”.
8. Click “Next”.
9. Click “Install”.  
The set-up program will begin copying the files. CS-S20 and the instrument USB drivers are installed in this order. This may take several minutes. Please wait.
10. This completes the installation process. Click “Complete”.

## 1.2 Creating a shortcut

When CS-S20 has been installed, a shortcut to CS-S20 is registered in the Windows Start Menu (for Windows 8 and Windows 8.1, the shortcut is placed on the Start screen). To place a shortcut in another location, create it with Windows Explorer or another program.

## 1.3 To uninstall

### [Windows 7]

1. Click the “Start” button. Select [Settings - Control Panel] and double-click the [Add/Remove Programs] icon.
2. Select [CS-S20] from the list of [Change or Remove Programs] and click the “Change/Remove” button.
3. A dialog box that enables you to change, correct or delete programs opens. Select [Delete] and click “Next”.
4. When the file deletion confirmation dialog box opens, click “OK”. The uninstall program will start.
5. When the uninstall process is completed, click “Complete”.

### [Windows 8]

1. Right-click in the lower-left corner of the screen.
2. Click [Programs and Features] on the menu that appears.
3. Select [CS-S20], and then click “Uninstall”.
4. A dialog box for confirming that you want to uninstall the application appears. Click “Yes”.
5. A dialog box for confirming that you want to update the software appears. Follow the instructions on the screen to complete the uninstall process.

### [Windows 8.1 and Windows 10]

1. Right-click on the Start button in the lower-left corner of the screen.
2. Click [Programs and Features] on the menu that appears.
3. Select [CS-S20], and then click “Uninstall”.
4. A dialog box for confirming that you want to uninstall the application appears. Click “Yes”.
5. A dialog box for confirming that you want to update the software appears. Follow the instructions on the screen to complete the uninstall process.

## 2. Connection

Connecting an instrument to your PC enables you to control the instrument with the CS-S20 software or to exchange data between the PC and instrument. Connection to the PC is performed with a USB cable.

After the instrument has been connected, follow the procedure in **2.2 Checking the COM port number** to determine which COM port number has been assigned to the instrument.

### Notes:

- Attach the connectors securely in the correct orientation.
- Do not touch the terminals of the connector. Keep them clean and do not apply excessive force.
- Use a cable of sufficient length. If the cable is too short, a connection failure or disconnection may result.
- To disconnect the cable, grip the plug or connector securely. Do not pull the cable or bend it with excessive force.

### 2.1 Connecting to a PC

#### 2.1.1 When the USB cable is used

1. Turn on both the instrument and the computer.
2. Connect the instrument and the computer via USB cable. Connect the computer end of the cable to the USB port of the computer. When the instrument is connected to the PC with a USB cable, and CS-S20 is activated, the remote communication mode will be automatically set.
3. When the instrument is connected to the computer for the first time, the USB driver will be automatically installed. Installation may take several minutes. When driver installation has been completed, the computer will automatically assign a COM port to the instrument.

### 2.2 Checking the COM port number

#### [Windows 7]

From the Start menu, right-click “Computer” and select “Properties” to open Device Manager. Click “Ports (COM & LPT)” to expand it, and the assigned COM port number is displayed.

#### [Windows 8]

Point to the lower-left corner of the screen, and then right-click in the window that appears. In the menu that appears, click “Device Manager” to open Device Manager. Click “Ports (COM & LPT)” to expand it, and the assigned COM port number is displayed.

#### [Windows 8.1 and Windows 10]

Point to the lower-left corner of the screen, and then right-click on the “Start” button that appears. In the menu that appears, click “Device Manager” to open Device Manager. Click “Ports (COM & LPT)” to expand it, and the assigned COM port number is displayed.

#### Note (All OS):

If the connected instrument is not shown in “Ports (COM & LPT)” and appears as “Unknown device” with an exclamation mark (“!”), continue from the above procedure by performing the procedure in **2.2.1 Manually updating the driver**.

#### 2.2.1 Manually updating the driver

If the instrument is not shown in “Ports (COM & LPT)” and appears as “Unknown device” with an exclamation mark (“!”), right-click on the “Unknown device”, select “Update driver” and select the folder where CS-S20 has been installed.

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