Color Data Software

SpectraMagic DX

- En** Installation Guide
- Fr** Guide d'installation
- De** Installationsanleitung
- It** Guida all'installazione
- Po** Instrukcja instalacji
- Pt** Guia de instalação
- Ru** Инструкция по установке
- Es** Guía de instalación
- Tr** Yükleme Rehberi
- Cs** 安裝指南
- Ct** 安裝指南
- JP** インストールガイド

KONICA MINOLTA

English

Français

Deutsch

Italiano

Polski

Português

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Español

Türk

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Software License Agreement

The terms of the license agreement of SpectraMagic DX software are provided in the Software License Agreement dialog box displayed on-screen during the installation process. This software can be installed only if you agree to all the terms of the agreement.

Safety Precautions

Before you use SpectraMagic DX software, we recommend that you thoroughly read this manual as well as the instruction manuals of your PC and the instrument.

Every effort has been made to ensure the accuracy of the content of this software. However, should you have any questions or comments, please contact the nearest KONICA MINOLTA authorized service facility.

Notes on Installation

- SpectraMagic DX application software is designed to be used with the Windows 10, or Windows 11 operating system. Note that no operating system is included with this software. One of these operating systems must be installed on the PC before this software can be installed.

Notes Regarding USB Devices (Flash Memory, Protection Key)

- When plugging the USB device into your computer, be sure that it is in the correct orientation. Do not forcibly plug it in.
- Do not touch the contacts of the USB flash drive.
- After using the USB device, return it to its case and store in a safe place.
- Avoid exposing the USB device to rapid temperature changes and condensation.
- Avoid leaving USB device in locations where it may be exposed to high temperatures from direct sunlight or heaters.
- Do not drop the USB device or subject it to strong impact.
- Keep the USB device away from water, alcohol, paint thinners, and other such substances.

Disposal Method

- Make sure that this product (USB device) is either disposed of or recycled correctly in accordance with local laws and regulations.

Notes on Instruction Manual

The Instruction Manual is installed in PDF form during the software installation. Versions of the Instruction Manual in other languages are also installed.

- If the software was purchased on a USB flash memory device, the instruction manuals can be read on that device in the \Manual folder.

To view the instruction manual while using the software, click on  in the upper right of the software window and select [Open SpectraMagic DX Manual] from the menu bar.

Formal designations of application software used in this manual

(Designation in this manual)	(Formal designation)
Windows, Windows 10	Microsoft® Windows® 10 Pro Operating System
Windows, Windows 11	Microsoft® Windows® 11 Pro Operating System

Trademarks

- "Microsoft", "Windows", "Windows 10" and "Windows 11" are registered trademarks of Microsoft Corporation in the U.S.A. and other countries.
- "KONICA MINOLTA", the Konica Minolta logo and symbol mark, "Giving Shape to Ideas", and "SpectraMagic" are registered trademarks or trademarks of Konica Minolta, Inc.
- "Intel" and "Pentium" are registered trademarks of Intel Corporation in the U.S.A. and other countries.
- Other company names and product names mentioned in this manual are the registered trademarks or trademarks of their respective companies.

1. System Requirements

1.1 System Requirements

OS	Windows 10 Pro 32-bit Windows 10 Pro 64-bit Windows 11 Pro <ul style="list-style-type: none"> English, Japanese, German, French, Spanish, Italian, Portuguese, Russian, Turkish, Polish, Simplified Chinese, and Traditional Chinese versions of the above OS. The hardware of the computer system to be used must meet or exceed the greater of the recommended system requirements for the compatible OS being used or the following specifications.
Computer	PC equipped with a processor equivalent to Intel Core i5 2.7GHz or better
Memory	At least 2 GB (4 GB or more recommended)
Hard disk drive	20 GB of available hard disk space <ul style="list-style-type: none"> At least 10 GB of available disk space is required on the system drive (drive where the OS is installed) for database..
Display	Display hardware capable of displaying 1280 x 768 pixels / 16-bit color or better
USB port	Required for protection key if used. Not necessary for electronic license.
USB or serial port	Required for connection to instrument

1.2 Compatible Instruments

CM-M6, CM-25cG, CM-700d, CM-600d, CM-2600d, CM-2500d, CM-2500c, CM-2300d, CM-3700A, CM-3700A-U, CM-3700d, CM-3600A, CM-3610A, CM-5, CR-5

1.3 Language

Display language	English, Japanese, German, French, Spanish, Italian, Portuguese, Russian, Turkish, Polish, Simplified Chinese, and Traditional Chinese (Selectable after installation.)
-------------------------	---

2. Installing/uninstalling the software

SpectraMagic DX software is available on a USB flash memory or by downloading from the Konica Minolta website:

<https://www.konicaminolta.com/instruments/download/software/color/index.html>

- When downloading the software from the Konica Minolta website, be sure to download the SpectraMagic DX edition matching the Windows edition of your computer. The 32-bit edition can be installed only on computer systems with 32-bit editions of Windows, and the 64-bit edition can be installed only on computer systems with 64-bit editions of Windows.

Setup file name:

The name of the setup file is different depending on how the installation files were obtained.

- If the software was received on a USB flash memory, the name of the setup file will be "setup.exe" in the root directory of the USB flash memory.
- If the software was downloaded from the Konica Minolta website, the setup file will be contained in a *.zip file. In this case, extract all files to a folder on your computer before proceeding. The name of the setup file will be "setup_x86.exe" (for 32-bit Windows) or "setup_x64.exe" (for 64-bit Windows).

2.1 Installation procedure

- When installing the software, you must log on to the computer as a user with Administrator privileges.
 - If a previous version of SpectraMagic DX has been installed on the computer, uninstall the existing SpectraMagic DX version before installing the new version.
 - A dialog box for checking the User Account Control may be displayed multiple times during the process as the software and various drivers are installed. When this dialog box is displayed, click "Continue".
 - If there are any protection keys (including protection keys for other software) connected to the computer, unplug them before proceeding with installation.
 - If there are any Konica Minolta instruments connected to the computer, unplug them before proceeding with the installation.
1. Start Windows on your computer.
 - If Windows Update is running, wait until Windows Update has finished and then reboot the computer before installing SpectraMagic DX.
 2. Browse to the setup file.
 - » When installing from a USB flash memory, browse to the root folder of the USB flash memory.
 - » When installing from a downloaded file, browse to the folder to which the downloaded zip file was extracted.
 3. Double-click on the setup file. The setup will start.
 - It may take a few minutes for the setup dialog to appear while the installer checks the current system environment.
 4. After a few seconds, a dialog for selecting the language to use during installation will appear.
 - This is the language to use during installation only. The display language for SpectraMagic DX can be selected after installing SpectraMagic DX by selecting *Application Settings* from the **Document** menu and selecting the language in the Start options category of the Application

Settings dialog.

- To select a different language, click on the current language and select the desired language for installation from the pull-down list of languages that appears
5. Click [OK]. The installation will proceed.
 6. Depending on the computer environment and currently installed software, the SpectraMagic DX installer may need to install certain necessary Microsoft components, such as SQL, .NET 4.5.2 Framework, etc. If the installer determines that Microsoft components need to be installed, the installer for those components will start automatically.
 - Depending on the Microsoft component to be installed, it may be necessary to agree to the License Agreement for that component.
 - A command prompt window will appear during the installation of SQL. DO NOT CLOSE THIS WINDOW. When SQL installation has finished, the command prompt window will close automatically.
 - In some cases, it may be necessary to reboot your computer after installing Microsoft components. In such case, reboot your computer and then restart the SpectraMagic DX installation program from step 1 above.
 - Installation may require 30 minutes or longer.
 7. Click [Next]. The SpectraMagic DX license will be shown.
 - To print the license, click [Print]. The license will also be stored as a file in the SpectraMagic DX installation folder.
 8. Click on [I accept the terms of the license agreement] to select it and enable the [Next] button, and then click [Next].
 9. A dialog showing the folder where SpectraMagic DX will be installed will be shown. If the default folder is acceptable, click [Next].
 - To change the folder, click [Change ...] and browse to the desired folder before clicking [Next].
 10. Click [Install]. The SpectraMagic DX installation will proceed.
 - Installation may require several minutes. Please do not do anything else on the computer until SpectraMagic DX installation is completed.
 - A dialog for driver installation will appear during installation of SpectraMagic DX. Be sure to install the driver. If canceled, installation of SpectraMagic DX will also be canceled.
 11. When SpectraMagic DX installation has been completed, click [Finish] in the dialog that appears to finish the installation and close the dialog.

2.1.1 Creating a shortcut

When SpectraMagic DX has been installed, a shortcut to SpectraMagic DX is registered in the Windows Start Menu. To place a shortcut in another location such as the Desktop, create it with Windows Explorer or another program.

2.2 Uninstallation procedure

Uninstallation of SpectraMagic DX is performed using the standard Windows uninstallation method.

- The database used by SpectraMagic DX will not be uninstalled when SpectraMagic DX software is uninstalled.
 - Microsoft components (SQL, etc.) were installed during the installation of SpectraMagic DX will not be uninstalled when SpectraMagic DX software is uninstalled. Since the Microsoft components may be used by other software, they should not be uninstalled.
 - In particular, DO NOT UNINSTALL THE SQL FEATURE "SMDXEXPRESS". Uninstalling this SQL feature will cause the database containing SpectraMagic DX data, etc. to become unusable, resulting in the loss of data.
1. Right-click on the Start button in the lower-left corner of the screen.
 2. Click [Programs and Features] on the menu that appears.
 3. Select [SpectraMagic DX], and then click "Uninstall".
 4. A dialog box for confirming that you want to uninstall the application appears. Click "Yes".
 5. A dialog box for confirming that you want to update the software appears. Follow the instructions on the screen to complete the uninstall process.

3. Connection

For information on physical connections to instruments, refer to the Instruction Manual for each instrument.

For information on how to make serial port settings and connect to the instrument from within SpectraMagic DX, refer to SpectraMagic DX Instruction Manual.

4. About the SpectraMagic DX End User License

The end user license for SpectraMagic DX is available in either hardware or electronic form; further, the electronic license is available as a single license installed on a single computer, or as a multi-user floating license. Please see the section below corresponding to the license form that you purchased.

4.1 Hardware protection key license

If you purchased SpectraMagic DX with a hardware license, you will have received a USB hardware protection key with your software.

When using SpectraMagic DX, the USB protection key must be plugged into the USB port of the computer. The license is active only while the USB protection key is plugged into the computer.

4.2 Electronic license

If you purchased SpectraMagic DX with an electronic license, you will have received a License Certificate (either as a paper certificate or a pdf file) that lists your Product Key.

- The Product Key is necessary for performing license activation. Please keep it in a safe place.

The license will not be valid until you have activated the license.

- Be sure that e-mails from webmaster@selms-mail.konicaminolta.com are not blocked by your firewall and are not labeled as spam.

4.2.1 Online activation

To perform online activation of your license, follow the procedure below.

- The computer must be connected to the internet to perform online activation.

1. Start SpectraMagic DX.
2. When the SpectraMagic DX Product Activation dialog appears, click [Online Activation].
3. The SpectraMagic DX Product Activation dialog will change to show a textbox for inputting your Product Key.
4. Input your Product Key (from your license certificate) into the space provided and click [ACTIVATE].
5. The software will contact the SE LMS website and the SpectraMagic DX Product Activation dialog will show the message "SpectraMagic DX License activation is in Progress" with a progress bar.
6. A message requesting that user registration be performed will be shown. Click [Open]. The "User information and license confirmation" screen of the Konica Minolta Sensing Software License Management System (SE-LMS) website will be shown.
 - If desired, the language of any screen in the SE-LMS website can be changed by clicking on Select Language at the upper left of the screen and selecting the desired language from the pulldown that appears.
7. Check that the license information shown in the upper part of the screen matches your order. If it matches your order, click on "This matches to the items of the order" to put a check in the checkbox.
 - If the information does not match your order, contact the salesperson from whom you purchased the software.

8. Fill in your information in the lower part of the screen.
 - At a minimum, your e-mail address (the e-mail address you want to use for user registration), first name, last name, company name and country are required information.
9. If you agree with the site's privacy policy, click on the checkbox next to "I agree about the handling of personal information" to put a check in the checkbox.
 - To see the privacy policy for this website, click on "I agree about the handling of personal information". A separate window showing the privacy policy will open.
 - You must agree with the privacy policy in order to complete the registration process.
10. Click [Confirmation and registration]. A confirmation screen will appear.
11. If this is your first time to register, create a password by inputting it in the 2 textboxes. If you have registered previously, input your password in the textbox.
12. Click [Registration]. The user registration process is complete, and an email confirming your registration will be sent to you.
13. After a few seconds, activation will be completed and SpectraMagic DX will restart. License activation is complete.

4.2.2 Offline activation

Offline activation can be performed on a computer that does not have internet access provided that another computer that does have internet access is available.

To perform offline activation of your license, follow the procedure below.

1. Start SpectraMagic DX. .
2. When the SpectraMagic DX Product Activation dialog appears, click [Offline Activation]. The dialog will switch.
3. Click [Create C2V file]. A Select Folder dialog will appear.
4. Select the folder in which to create the C2V file and click OK. The file will be created.
5. Copy that file to a computer which has internet access.
6. Go to the SE LMS website at the following URL:
<https://selms.konicaminolta.com/login>
 - A link to the website is also provided on the License Information category screen of the Application Settings dialog of SpectraMagic DX.
7. The SE-LMS License authentication screen will be shown. Input your Product Key (from your license certificate) into the space provided and click [SEND].
8. Input your email address and password, and click [Log in]. The SE-LMS screen will be shown, with a list of licenses for your registered email address. To see the full list, click on the "+" at the upper left of the list.
9. In the list of licenses, locate the Product Key for which you want to perform offline activation and click [Offline Activation] for that product key.
10. A screen showing the selected Product Key and requesting the C2V file created in step 4 will appear. Click [Choose File] to open the Open dialog, browse to the C2V file and click [Execute].
 - If you wish to have the resulting V2C file sent to a different e-mail address, input that e-mail address in the space provided before clicking [Execute].
11. An e-mail confirming completion of activation will be sent to your e-mail address (or alternate

specified e-mail address) and the screen will change to Complete the Off-line activation, with a [Download V2C file] button.

- 12.** Click [Download V2C file] and save the file on your computer.
- 13.** Copy the downloaded V2C file to the computer for which the license will be activated.
- 14.** in the SpectraMagic DX Product Activation dialog (as shown in step 2), click [Browse V2C file] and browse to the copied V2C file.
- 15.** Click [ACTIVATE].
- 16.** The SpectraMagic DX Product Activation dialog will show the message "SpectraMagic DX License activation is in Progress" with a progress bar.
- 17.** After a few seconds, activation will be completed and SpectraMagic DX will restart. License activation is complete.

4.2.3 Using a Floating License

The floating license is a multi-user license that is installed on a single computer on a network (which acts as the license server) and can then be used by other computers (client users) on the network up to the number of licenses purchased.

- The license server PC and the client user PCs must be on the same subnet. (If you wish to use a floating license with PCs on a different subnet, please contact your Konica Minolta sales representative.)
- The license server PC must be on while a client PC is using a license.
- For details on the range of your subnet, please check with your IT administrator.

To install, activate, and use a floating license, please follow the procedure below.

1. Install SpectraMagic DX on the PC that will be used as the license server.
2. Start SpectraMagic DX.
3. Activate the license installed on the license server PC according to the procedure in "4.2.1 Online activation" or "4.2.2 Offline activation".
4. Install SpectraMagic DX on the PCs which will be client PCs.

The floating license is now activated. When SpectraMagic DX is started on a client PC, it will connect to the license server PC to obtain a license. The number of PCs which can use SpectraMagic DX at the same time will be the number of user licenses in the floating license.

Restricting license connection range

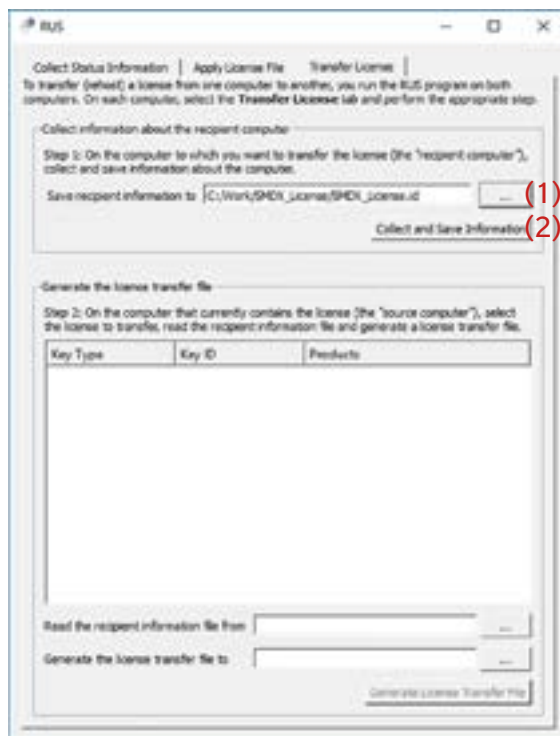
The license connection range of the license server and client PC can be performed by connecting to <http://localhost:1947> from each PC and using the Admin Control Center. For details on how to use the Admin Control Center, refer to the Admin Control Center help.

- Default setting: All PCs which are on the same subnet as the license server PC and which have SpectraMagic DX installed are within the connection range.

4.2.4 License transfer

To transfer (rehost) a license from one computer to another, you run the RUS program on both computers. On each computer, select the **Transfer License** tab and perform the procedure below.

- In advance, SpectraMagic DX must be installed on the recipient computer (the computer to which the license is transferred).
1. Open the RUS tool on the recipient computer.
 - Path: C:\Program Files\KONICA MINOLTA\SpectraMagic DX
 - Name: RUS_100634.exe
 2. Create a recipient file (id file) on the recipient computer.
 - (1) Specify a path and filename you like.
 - (2) Click this button. Then id file is created.
 3. Copy the id file from the recipient computer to the source computer (the computer from which the license will be transferred).

4. Open the RUS tool on the source computer.
5. Create a license transfer file (h2h file) on the source computer.
 - (3) Specify the id file copied from the recipient computer before.
 - (4) Specify a path and filename of h2h file you like.
 - (5) Click this button. Then h2h file is created.

6. Copy the h2h file to the recipient computer.

7. Apply a license transfer file (h2h file) on the recipient computer, and the license will be transferred.

- (6) Change to this tab.
- (7) Specify the created h2h file.
- (8) Click this button. Activation is complete.

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