# Color Data Software

# SpectraMagic NX2

- En** Installation Guide
- Fr** Guide d'installation
- De** Installationsanleitung
- It** Guida all'installazione
- Po** Instrukcja instalacji
- Pt** Guia de instalação
- Ru** Инструкция по установке
- Es** Guía de instalación
- Tr** Yükleme Rehberi
- Cs** 安裝指南
- Ct** 安裝指南
- Jp** インストールガイド
- Ko** 설치 가이드

KONICA MINOLTA

English

Français

Deutsch

Italiano

Polski

Português

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Español

Türk

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## Software License Agreement

The terms of the license agreement of SpectraMagic NX2 software are provided in the Software License Agreement dialog box displayed on-screen during the installation process. This software can be installed only if you agree to all the terms of the agreement.

### Safety Precautions

Before you use SpectraMagic NX2 software, we recommend that you thoroughly read this manual as well as the instruction manuals of your PC and the instrument.

Every effort has been made to ensure the accuracy of the content of this software. However, should you have any questions or comments, please contact the nearest KONICA MINOLTA authorized service facility.

## Notes on Installation

- SpectraMagic NX2 application software is designed to be used with the Windows 10, or Windows 11 operating system. Note that no operating system is included with this software. One of these operating systems must be installed on the PC before this software can be installed.

## Notes Regarding USB Devices (Flash Memory, Dongle)

- When plugging the USB device into your computer, be sure that it is in the correct orientation. Do not forcibly plug it in.
- Do not touch the contacts of the USB flash drive.
- After using the USB device, return it to its case and store in a safe place.
- Avoid exposing the USB device to rapid temperature changes and condensation.
- Avoid leaving USB device in locations where it may be exposed to high temperatures from direct sunlight or heaters.
- Do not drop the USB device or subject it to strong impact.
- Keep the USB device away from water, alcohol, paint thinners, and other such substances.

## Disposal Method

- Make sure that this product (USB device) is either disposed of or recycled correctly in accordance with local laws and regulations.

## Notes on Instruction Manual

The Instruction Manual is installed in PDF form during the software installation. Versions of the Instruction Manual in other languages are also installed.

- If the software was purchased on a USB flash memory device, the instruction manuals can be read on that device in the \Manual folder.

To view the instruction manual while using the software, press the F1 key or select Help - Manual ...

## Formal designations of application software used in this manual

(Designation in this manual)	(Formal designation)
Windows, Windows 10	Microsoft® Windows® 10 Pro Operating System
Windows, Windows 11	Microsoft® Windows® 11 Pro Operating System

## Trademarks

- "Microsoft", "Windows", "Windows 10" and "Windows 11" are registered trademarks of Microsoft Corporation in the U.S.A. and other countries.
- "KONICA MINOLTA", the Konica Minolta logo and symbol mark, "Giving Shape to Ideas", and "SpectraMagic" are registered trademarks or trademarks of Konica Minolta, Inc.
- "Intel" and "Pentium" are registered trademarks of Intel Corporation in the U.S.A. and other countries.
- Other company names and product names mentioned in this manual are the registered trademarks or trademarks of their respective companies.

# 1. System Requirements

---

## 1.1 System Requirements

<b>OS</b>	Windows 10 Pro 64-bit Ver. 1903 or higher Windows 11 Pro <ul style="list-style-type: none"> <li>English, Japanese, German, French, Spanish, Italian, Portuguese, Russian, Turkish, Polish, Simplified Chinese, Traditional Chinese, and Korean versions of the above OS.</li> <li>The hardware of the computer system to be used must meet or exceed the greater of the recommended system requirements for the compatible OS being used or the following specifications.</li> </ul>
<b>Computer</b>	PC equipped with a processor equivalent to Intel Core i5 2.7GHz or better
<b>Memory</b>	2 GB (4 GB or more recommended)
<b>Hard disk drive</b>	10 GB of available hard disk space
<b>Display</b>	Display hardware capable of displaying 1280 x 768 pixels / 16-bit color or better
<b>USB port</b>	Required for dongle version. Not necessary for electronic license.
<b>Connection to external network</b>	Required for license activation

## 1.2 Compatible Instruments

CM-36dGV, CM-36dG, CM-36d, CM-26dG, CM-26d, CM-25d, CM-23d, CM-25cG, CM-3700A, CM-3700A-U, CM-3600A, CM-3610A, CM-M6, CF-300, CM-5, CR-5, CM-17d, CM-16d, CM-512m3A, CM-700d, CM-700d-U, CM-600d, CM-2600d, CM-2500d, CM-2500c, CR-400, CR-410, DP-400

## 1.3 Language

<b>Display language</b>	English, Japanese, German, French, Spanish, Italian, Portuguese, Russian, Turkish, Polish, Simplified Chinese, Traditional Chinese, and Korean (Selectable after installation.)
-------------------------	---

## 2. Installing/uninstalling the software

---

SpectraMagic NX2 software is available on a USB flash memory or by downloading from the Konica Minolta website:

<https://www.konicaminolta.com/instruments/download/software/color/index.html>

### Setup file name:

The name of the setup file is different depending on how the installation files were obtained.

- If the software was received on a USB flash memory, the name of the setup file will be "setup.exe" in the root directory of the USB flash memory.
- If the software was downloaded from the Konica Minolta website, the setup file will be contained in a \*.zip file. In this case, extract all files to a folder on your computer before proceeding.

### Configuration Tool CM-CT1:

The setup file does not include Configuration Tool CM-CT1. Configuration Tool CM-CT1 can be downloaded from the Konica Minolta website:

<https://www.konicaminolta.com/instruments/download/software/color/index.html>

## 2.1 Installation procedure

- When installing the software, you must log on to the computer as a user with Administrator privileges.
  - If a previous version of SpectraMagic NX2 has been installed on the computer, uninstall the existing SpectraMagic NX2 version before installing the new version.
  - A dialog box for checking the User Account Control may be displayed multiple times during the process as the software and various drivers are installed. When this dialog box is displayed, click "Yes".
  - If there are any dongles (including dongles for other software) connected to the computer, unplug them before proceeding with installation.
  - If there are any Konica Minolta instruments connected to the computer, unplug them before proceeding with the installation.
1. Start Windows on your computer.
    - If Windows Update is running, wait until Windows Update has finished and then reboot the computer before installing SpectraMagic NX2.
  2. Browse to the setup file.
    - » When installing from a USB flash memory, browse to the root folder of the USB flash memory.
    - » When installing from a downloaded file, browse to the folder to which the downloaded zip file was extracted.
  3. Double-click on the setup file. The setup will start.
    - It may take a few minutes for the setup dialog to appear while the installer checks the current system environment.
  4. After a few seconds, a dialog for selecting the language to use during installation will appear.
    - This is the language to use during installation only. The display language for SpectraMagic NX2 can be selected after installing SpectraMagic NX2 by opening the Application Settings of the SpectraMagic NX2 Launcher and selecting Display Language in the User Settings category.

- To select a different language, click on the current language and select the desired language for installation from the pull-down list of languages that appears
- 5.** Click [ OK ]. The installation will proceed.
  - 6.** Depending on the computer environment and currently installed software, the SpectraMagic NX2 installer may need to install certain necessary Microsoft components, such as NET 4.5.2 Framework, etc. If the installer determines that Microsoft components need to be installed, the installer for those components will start automatically.
    - Depending on the Microsoft component to be installed, it may be necessary to agree to the License Agreement for that component.
    - In some cases, it may be necessary to reboot your computer after installing Microsoft components. In such case, reboot your computer and then restart the SpectraMagic NX2 installation program from step 1 above.
    - Installation may require 30 minutes or longer.
  - 7.** Click [ Next ]. The SpectraMagic NX2 license will be shown.
    - To print the license, click [ Print ]. The license will also be stored as a file in the SpectraMagic NX2 installation folder.
  - 8.** Click on [ I accept the terms of the license agreement ] to select it and enable the [ Next ] button, and then click [ Next ].
  - 9.** A dialog showing the folder where SpectraMagic NX2 will be installed will be shown. If the default folder is acceptable, click [ Next ].
    - To change the folder, click [ Change ... ] and browse to the desired folder before clicking [ Next ].
  - 10.** Click [ Install ]. The SpectraMagic NX2 installation will proceed.
    - Installation may require several minutes. Please do not do anything else on the computer until SpectraMagic NX2 installation is completed.
    - A dialog for driver installation will appear during installation of SpectraMagic NX2. Be sure to install the driver. If canceled, installation of SpectraMagic NX2 will also be canceled.
  - 11.** When SpectraMagic NX2 installation has been completed, click [ Finish ] in the dialog that appears to finish the installation and close the dialog.
  - 12.** If you have a SpectraMagic NX2 license, perform license activation according to the License Activation procedure in "1.1.3.1 License Info" of the SpectraMagic NX2 Instruction Manual.
    - If you do not have a license and this is the first time SpectraMagic NX2 has been installed on this computer, you will be able to use all functions for a trial period of 30 days. After the trial period has ended, it will be necessary to purchase a license to continue using the software.

## 2.2 Uninstallation procedure

Uninstallation of SpectraMagic NX2 is performed using the standard Windows uninstallation method.

- Microsoft components were installed during the installation of SpectraMagic NX2 will not be uninstalled when SpectraMagic NX2 software is uninstalled. Since the Microsoft components may be used by other software, they should not be uninstalled.
1. Right-click on the Start button in the lower-left corner of the screen.
  2. Click [Programs and Features] on the menu that appears.
  3. Select [SpectraMagic NX2], and then click "Uninstall".
  4. A dialog box for confirming that you want to uninstall the application appears. Click "Yes".
  5. A dialog box for confirming that you want to update the software appears. Follow the instructions on the screen to complete the uninstall process.

## 3. Connection

---

For information on physical connections to instruments, refer to the Instruction Manual for each instrument.

For information on how to connect to the instrument from within SpectraMagic NX2, refer to SpectraMagic NX2 Instruction Manual.

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